Love of life begins with "Rcom"

Ver. 1.0

USER'S MANUAL

INSECT MAX 70






Rcom Insect Incubator Max 70 User Manual

Contents





How to Use Rcom Insect MAX 70

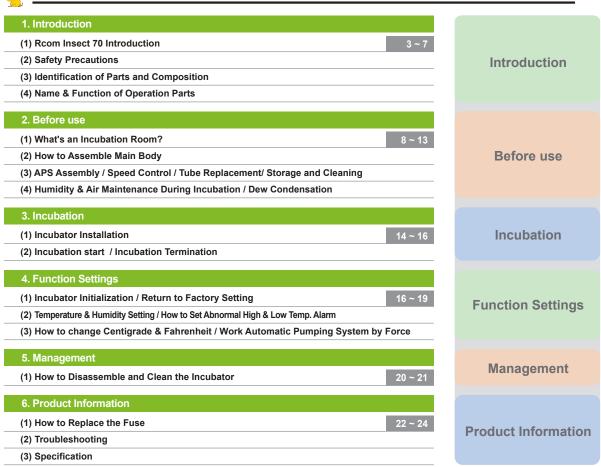

<u>^</u>

Successful hatching depends on many numbers of factors. Neither Autoelex (Rcom) nor its global distributors and vendors can be held responsible for loss of life or property damage caused by user neglect, alteration, modification or painting, change of use or power failure. To avoid errors, mishaps and danger, carefully check this User Manual before initial use. Also please check the manual regularly to ensure your operation of the machine is correct.



2



1. Introduction

(1) Rcom Insect 70 Introduction

Rcom Insect 70

First of all, we would like to express our deep appreciation for all the customers who purchased Rcom Insect Incubator. Rcom Insect Incubator was designed to be used easily and conveniently by the newest digital control method for controlling optimal incubation environment automatically, however, in order to get the best incubation result, at first, it is very important for users to have some incubation knowledge and right machine operation method. Therefore, before suing incubator, be sure to read instruction manual. Rcom Insect Incubator provides the best incubation environment, however, it is necessary for users to observe and adjust each environment necessary for incubation including egg kind, temperature and humidity management by incubation date, surrounding conditions of incubator, and this product was made it possible to make the best setting for all the environments necessary for incubation.

Features of Rcom Insect 70

[Main Features]

- * Luxury design with streamlined egg shaped concept
- * Automatic Temperature&Humidity setting and control
- * Artificial intelligence electronic control device to automatically adjust the cycle of incubator according to the ambient conditions
- * Convenient humidifier with Automatic Pumping System (A.P.S)
- * Clear and transparent large view-window
- * Minimizing the parasitic bacteria by not using any interior material and using revolutionary hygienicduplicated insulation structure (bottom tray)
- * Variable air control lever to control air amount inside the incubator
- * Applying Rcom's optimum air flow technology to avoid the fan's air from directly touching egg
- * Enhanced reliability by applicating Swiss's Sensirion's 3rd generation temperature & humidity sensor

[Easy Features]

- * Degree C / degree F interchangeable
- * Alarm and indicator function for abnormal high or low temperature by drastic and unusual outer temperature change
- * Incubation data memory function and power outage notification function in case of power outage
- * Closed structure to avoid waterdrops (condensation) on the view window from leaking out of incubator
- * Rotating heater support which conveniently controls the tension of the heater
- * Application of Water Nipple to supply water easily for humidification
- * Automatic humidification function for minimum 2 minutes (Press + button for 10 sec.)
- * Automatic Humidification Pumping System, and precise Speed Regulator VR to control the amount of pumping
- * Air Vent Holes for inflow of fresh air from outside to inside by affecting the heat insulation to the minimum
- * Not required to have a balance weight with the optimized incubator design considering the centroid of incubator's top and bottom when eggs are placed
- * The Automatic Pumping System (APS) for accurate humidification, as a pump of silicone tube, is structured to make tube changed with easy and durable by installing mini roller on the four areas with friction

1. Introduction

(2) Safety Precautions

INSECT TO

Electrical hazards

ACAUTION Be careful the details below when you use.



Do not use a damaged power cord or loose outlet.

Risk of electric shock or fire.



Do not pull the cord when taking out the power cord, and outlet during the incubation keep wet hands away from connecting plug.

▶ Risk of electric shock or fire



Never pull the plug out of the

Incubation will be interrupted



Do not twist or crush electric cord.

Risk of electric shock or fire.



Do not insert multiple connecting plugs in an outlet.

Risk of fire or electrical

Setting Cautions

ACAUTION Be careful the details below when you use.



Do not install in dusty or dirty environment

Risk of damage or fire within the incubator



Do not install under the direct sunlight.

▶ Risk of fire or interference with the incubation process



Do not install in moist or humid environment

Risk of fire or electric shock



Do not install in excessively cold or hot conditions. cigarette smoke, etc.

Risk of interference with the incubation process



Do not use any other non-standard parts except those provided.

Risk of damage or hatching failure.



Do not cover the ventilating opening.

► Inner temperature can rise, interrupting incubation



Install away from heat sources.

▶ Risk of damage to the incubator case and interference with the incubation process



Ensure that the incubator is installed on a stable surface away from edges.

 Risk of damage to incubator and eggs and user from accidental knocks or drops



Do not turn the incubator upside down.

Water wil pour out of the incubator, and view window can be dropped or damaged.



Children should be supervised to ensure that they do not play with the appliance.

Risk of knocking the incubator or accidental interference with the controls.



Do not disassemble or modify the incubator in any way.

Risk if electric shock or fire.



Please ensure that no small objects get into the holes on the incubator.



If the incubator sounds strange or emits smoke contact your service center.



Clean the incubator thoroughly before storing



If the incubator requires repair disconnect from the power supply and please contact to your service center.

Cleaning Cautions

ACAUTION

Be sure to disconnect the electric cord from the outlet before cleaning.



on the incubator surface.



cloth with a neutral detergent.



Brush away dust on the plug with a dry cloth.



Do not use chemicals like wax, For special cleaning of inner benzene, alcohol, thinner, aromatic, or lubricant, etc



part per year, contact to the service center.

If you don't clean the inside of


incubator for a long time, dust can cause some trouble or a fire.

INSECT TO

1. Introduction

(3) Identification of Parts and Composition


Identification of Parts

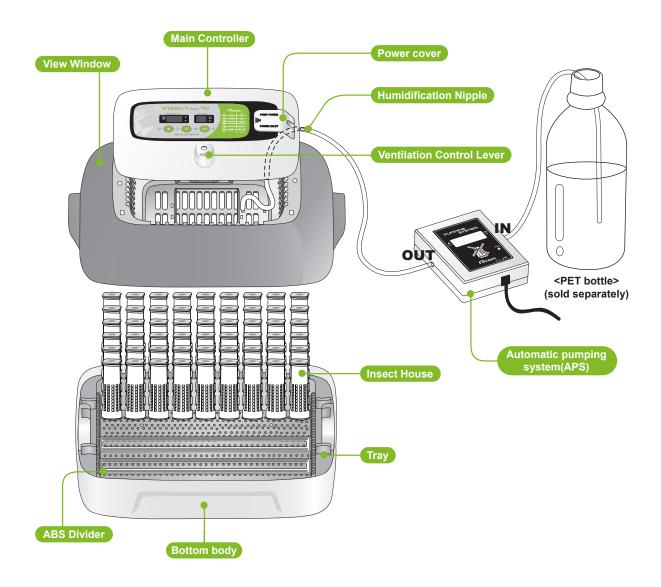

1. Introduction

(3) Identification of Parts and Composition



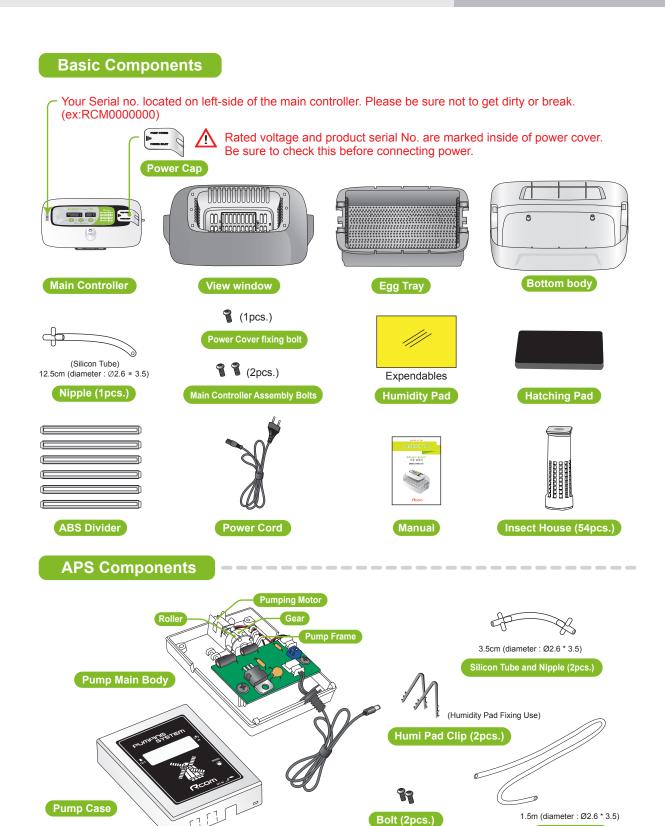

6

DIGITAL INCUBATOR


Silicon Tube

INSECT TO

1. Introduction

(4) Name & Function of Operation Parts

Name & Function of Operation Parts

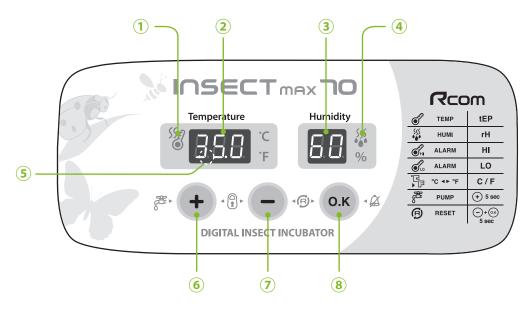

Names and components of each part

- ① Heater Operation Lamp
 - : Light on when working
- 2 Temp. Display
 - : Current Temperature Display
- ③ Humidity Display
 - : Current Humidity Display
- **4** Pumping System Operation Lamp
 - : Light on when working

- **5** Power Failure Alarm Function (FND light)
 - : Light on when power failure (Cancel Button : OK)
- 6 Up Button
 - : Setting Temp. & Humidity UP
- **7** Down Button
 - : Setting Temp. & Humidity DOWN
- **8 Menu Selecting Button**
 - : Setting Value check
- Insect MAX 70 is available to automatic humidity raise only, if you want to control low humidity level, please adjust "Air Controlling Lever" by manual [refer to page 13]

+ ·8· -	Setting Mode: Press two buttons at the same time to go to Setting Mode.
<u>О.К</u>	Select Menu / Quick Movement / Cancellation / Setting value check during incubation
(+)	Value + / APS is forced to operate if press 5sec. / APS is forced to operate for 2min. if press 10sec.
<u> </u>	Value - / Factory Initialization

• Function Key Press (+) (-) buttons at the same time to go to setting mode.

Function	Temp.		Humidity		Abnormal high temp. alarm		Abnormal low temp. alarm		Centigrade / Fahrenheit
Dispaly	tEP	→	rH	-	HI	-	LO	-	°C / °F
Default Setting	35.0℃		60%		2°C		-3°C		°C

2. Before use

(1) What is an incubation Room? (2) How to Assemble

INSECT TO


What is an Incubation Room?

Incubation room? It means the place where the incubator is installed and operated, and as it has much effect on incubation rate according to the environment of incubation room, it is good to install incubator at the place where there is little noise and vibration in the vicinity and the temperature range isn't severe in the temperature of $22\sim25^{\circ}$ C (71.6~77°F). In particular, as there are many cases that the temperature falls rapidly in the nighttime differently from the daytime, it is necessary to check certainly, and during daytime, incubator must not be exposed directly to the sunlight.




When the temperature of incubation room is lower than, recomended setting temperature there could be water-droplets in side of incubation room or water leak on the floor. because of dew condensation.

How to Assemble

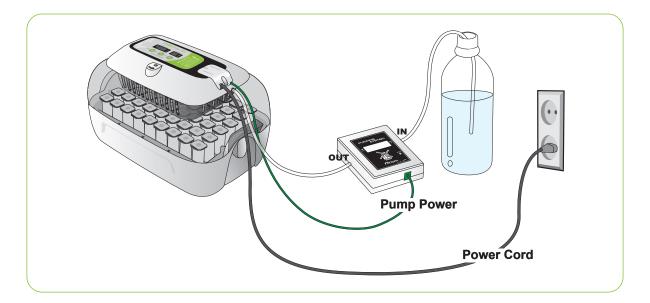

[How to Assemble Main Body]

- ▶ Please remove all packing from the incubator and parts. Then, check each part is present and undamaged. (Please keep the carton and packing materials for repacking to retain the incubator in good condition when not using.)
- ► (EX:RCM0000000) If you register your information with your machine's serial number on our web-site(www. Rcom.co.kr), you can get 2 year warranty service. [Refer to the page 24]

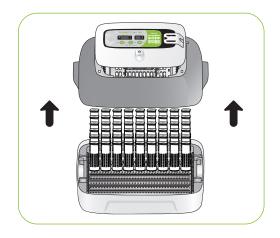


8

INSECT TO


2. Before use

(2) How to Assemble Main Body


▶ Loose two bolts connecting main controller and view window. After that, disassemble main controller from view window by pulling 4 hooks on main controller.

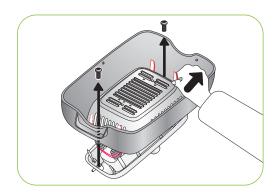

▶ Fix silicone tube which is assembled with nipple into the

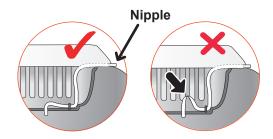


A Be careful to ensure the silicone tube is not folded.

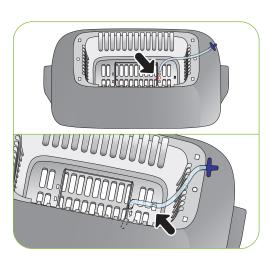

▶ Insert the nipple which is connected with view window to the hole of main controller, and assemble view window and main controller with two bolts.



Do not tighten the bolts too much.

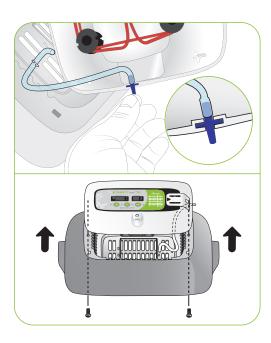



2. Before use

(2) How to Assemble Main Body

INSECT TO

Cut Humidity pad to proper size, and then fix on the view window with 2 pins included.

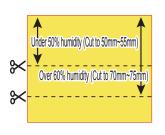

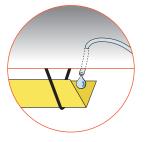

New Evaporating ability is different from pad size.

- You can use Humidity pad for about 4 ~ 6 months, but it can be different depending on water quality. (DISTILLED WATER is recommended to use.)
- * Humidity pad can be purchased separately.
- ▶ Assemble bottom body, Tray, and Tray hatching pad.





- ▶ Place Insect house by adjusting tray divider.
- ▶ Please close the view window.





Test the incubator to be sure that the incubator works without any problem before placing eggs in the incubator. Please ensure you remember how to use the machine and maybe re-read instructions.

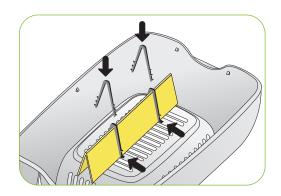

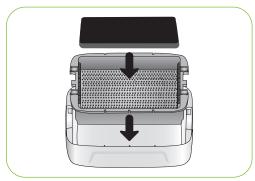






10

INSECT 70

2. Before use

(3) APS Assembly / Speed Control / Tube Replacement/ Storage and Cleaning

How to Assemble APS (Automatic Pumping System)

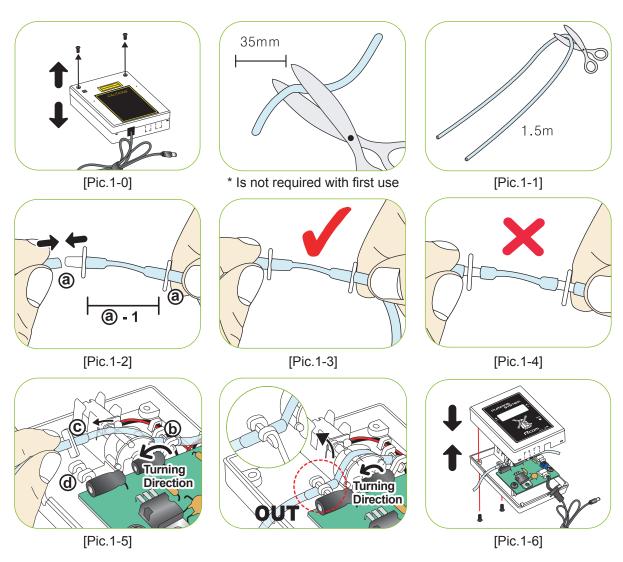

- ① Cut the silicone tube exactly to 35mm, and insert to nipple as picture 1-2. (When you purchase, product is originally assembled as @-1.)
- ② Cut the silicone tube of 1.5m to half, and insert to assembled nipple ③.
- ③ It is normal if you assembled as pic.1-3. If you do not insert silicon tube to the nipple completely as pic.1-4, pumping will not work well.
- (4) Loose two bolts of APS case as picture 1-0, and insert the assembled silicone tube and nipple as pic.1-3 to the side of pic. 1-5 (b).
- ⑤ Grip the part ⓒ on Pic.1-5, and pull the silicone tube so that fit to ⓓ. (If silicon tube should be assembled tightly.)
- ⑤ Spread IN and OUT of tube so that the silicone tube is not tangled. Insert tube on the groove, and cover the case. Here, be sure to check the silicone tube or wire is not jammed. Refer to the Pic.1-6.


2. Before use

(3) APS Assembly / Speed Control / Tube Replacement/ Storage and Cleaning



How to Control Water Flow Speed in APS

- The amount of water for humidification can be easily controlled with APS.
- With cross type screw driver, turn "Speed" clockwise to increase water flow and vice versa to decrease.
- ▶ To prevent water over flowing and dropping down in to the chamber, it is recommended to use the APS with lowest water flow speed. (1 drop of water in 3 sec.)

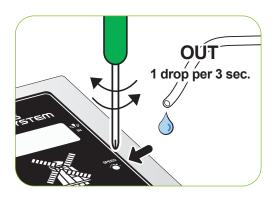



Humidity will raise up very slowly as artificial intelligence function check and memorize ambient conditions to keep incubator inside with optimum conditions.

How to Replace the Tube

▶ If the silicone tube wear out and it does not pump water effectively, you should replace silicone tube. [Refer to the page 13]



Be sure to use the silicon tube supplied from Autoelex(Rcom). (\emptyset 2.6 × 3.5)



If water does not flow out even if pumping motor is working, check if silicon tube is blocked or folded.

* If it failed in pumping water, check direction of IN and OUT again, and also check if motor is turning with anticlockwise. Also, length of silicon tube should be 35mm and inserted correctly. [Refer to the page 13]

Storage and Cleaning

- ▶ Do not wash APS with water but wipe with soft cloth.
- ▶ In case of long time storage, detach the silicon tube from the APS and remove the water from the tube. Keep the tube unkinked. It stops the tube hole from blocking.
- ▶ When assembling again, stretch the tube to open the hole in case the tube is blocked.

12

INSECT 70

2. Before use

(4) Humidity & Air Maintenance **During Incubation / Dew Condensation**

Humidity & Air Maintenance During Incubation

- ▶ When hatching, it's recommended not to open the lid often. If you open the lid often humidity will be rapidly decreased and it will take a long time to regain the proper humidity.
- * Air Controlling Lever : Outer fresh air can be flowed into incubator inside without affecting insulation. When eggs start hatching, open air controlling lever either fully or half-open.





 $\eta_{ip}^{(r)}$ Check water level every three days during incubation, and refill when needed. You'd better to use tap water than purified water or underground water for avoiding evaporating pad damage. (Evaporating pad is available from Autoelex Co.,Ltd. or Rcom distributors.)


Evaporating pad is expendables.

Dew Condensation

▶ This is a naturally occurring phenomenon when there is a significant difference in temperature between inside and outside of the incubator during incubation period. If this occurs, water may form in inside bottom of the incubator(bottom part).



No not place any goods near the incubator.

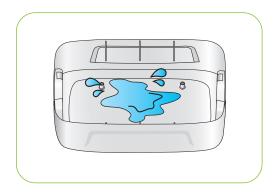

13

3. Incubation

(1) Incubator Installation



Incubator Installation


Test the incubator to be sure that the incubator works without any problem before placing eggs in the incubator. Please ensure you remember how to use the machine and maybe re-read instructions.

Open the power cap, and insert the incubator power cord and pump power cord.

Be careful to ensure the cords do not get jammed when you tighten screw bolt.




If you open power cap, there is a sticker which is showing rated current in the position of pic. a.

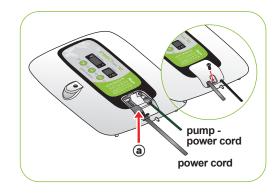

- ► Connect silicon tube end of APS (Automatic Pumping System) into the nipple of incubator, and the other end into PET bottle.
- ▶ Be sure to check rated current Pic. ⓐ before connecting power cord to outlet.
- ▶ Press (+) button for about 10 sec. for pump operation, and it automatically stop after 2min. If you want to stop pump operation, press any key.

If you do not insert silicon tube exactly, it will not work properly. [Refer to the page 11. How to replace silicon tube.]

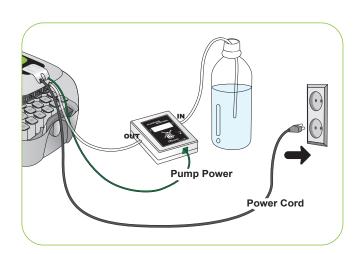

14



3. Incubation


(2) Incubation Start / **Incubation Termination**


How to Start Incubation


- If you connect power cord, incubator starts incubation with factory set conditions. [Factory setting : Temperature 35.0°C (95°F), Humidity RH 60%]
- Start incubation with required temperature & humidity for your species of birds.




There can be some odor the first time you use incubator, which is normal.

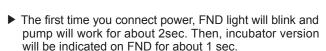

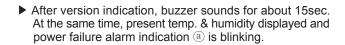


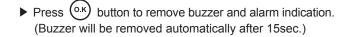

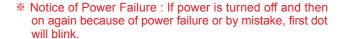



 $\eta_{ip}^{(i)}$ Quick Start : If you just connect power, it automatically starts incubation with factory setting. [Factory setting : Temperature 35.0°C (95°F), Humidity RH 60%]

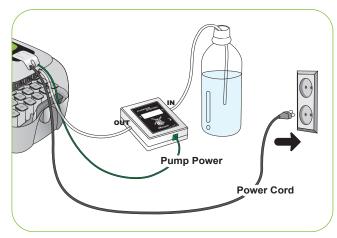





- Incubator display current temperature and humidity, and it goes to setting conditions within an hour.
- * The artificial intelligence system memorizes and classifies ambient conditions for keeping optimum temperature, and temperature can move up slowly at the first stage.

15

3. Incubation


4. Function Settings

(2) Incubation Start / Incubation Termination (1) Incubator Initialization / Return to Factory Setting




Incubation Termination

- ▶ There is no specified ON or OFF button on the incubator. Just turn off the power cord after hatching.
- It is not necessary to do Incubator Initialization with Insect MAX 70 incubator after hatching. Just connect power again when you want to start incubation.

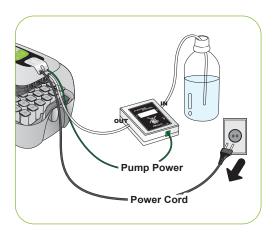

Incubator Initialization

- This function can be used when user input wrong setting, or incubation termination.
- ▶ Press (-) and (o.k) button at the same time for about 5sec. then display will show "rSt" and incubator returns to default setting.

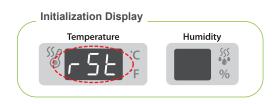




Calibrated setting value will not be initialized with this function. [If you need to initialize calibrated setting, refer to FACTORY SETTING next page.]

Return to Factory Setting

- ▶ This function is for return incubator to factory setting.
- ▶ Unplug power cord. Replug during button is press ed. Then, "rSt AL" will indicated in display, and incubator return to factory setting. [Default setting]



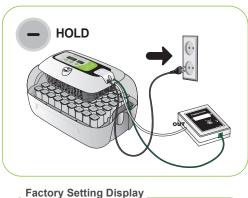



16



4. Function Settings

(2) Temperature Setting&Humidity Setting / How to Set Abnormal High Temp. Alarm

Temperature Setting




Function Key

Function	Temp.	Humidity	Abnormal high temp. alarm	Abnormal low temp. alarm	Centigrade / Fahrenheit
Dispaly	tEP =	rH •	→ HI •	→ LO •	→ °C / °F
Default Setting	35.0℃	60%	2°C	-3°C	°C

- ► Then, you can adjust temperature with


 and button. [Default setting: 35°C (95°F)]
- ▶ After setting required temperature, press ok button once then setting will be stored.

 Next, "rH" will indicated in humidity display for about 0.5 sec. and then humidity display will blink.
- Just press (ок) button if you do not need to change setting.

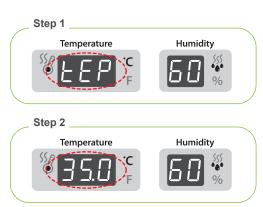

Humidity Setting

- ► From the temperature setting, press (o.k) button once and "rH" (Humidity) will be displayed for 0.5 second and humidity display will blink.
- ▶ When humidity display is blinking, adjust humidity setting with (+) or (-) button. [Default setting : RH 45%]
- ► After setting required humidity, press (o.k) button once then setting will be stored. Next, "HI" will be indicated for about 0.5sec.

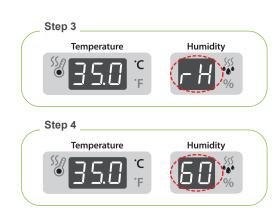



Insect MAX 70 is available to automatic humidity raise only, if you want to control low humidity level, please adjust "Air Controlling Lever" by manual [refer to page 13]

The temperature and humidity level is optimized according to the international standard at the time of ship ping. It could be recalibrated by the user if the user wishes to do so, but this is not recommended. When in need of adjustment, please refer to our homepage (www.Rcom.co.kr o Information o User Manual Category) or contact the place of purchase. Commercially available thermometer and hygrometers may have drastic measurement deviations, so it is recommended to use thermometer and hygrometer designated for this purpose. [Designated Thermometer and hygrometer – Refer to Rcom Homepage]

4. Function Settings

(2) Temperature Setting&Humidity Setting / How to Set Abnormal High Temp. Alarm



How to Set Abnormal High Temperature Alarm

* Max. abnormal temp. alarm limit (0.0 ~ 5.0°C Default Setting:2.0°C)



Abnormal High temperature:

This function gives notice when incubator temperature is higher than setting temperature because of abnor

mal ambient temperature increase.

When room temperature is higher than incubator tem perature, this function works with buzzer. Here, "HI" and temp. gap are displayed alternatively. Press OK button to remove buzzer.

- ► From the humidity setting, press the (o.k) button once and "HI" (abnormal HIgh temperature alarm) will be displayed for 0.5 second and temperature display will blink.
- ► After setting required value, press (o.k) button to save. Then, it will go to the next stage, abnormal low temp. setting with display "LO" for about 0.5 sec.





ex:





- It shows that the set temperature value is approximately 3°C (3°F) higher than the Maru incubator's incubation chamber(the Cabinet).
- * Just press (o.k) button if you do not need to change setting.

How to Set Abnormal Low Temperature Alarm

- ➤ This function gives notice when incubator temperature is lower than setting temperature because of abnormal ambient temperature decrease.
- From the Abnormal High Temperature Alarm setting, press the ox button once and "LO" (abnormal LOw temp. alarm) will be displayed for 0.5 second and the temperature display will blink.
- ► Adjust abnormal low temp. setting with to or button. [Default Value: -3°C (-3°F)]
- ► After setting required value, press OK button to save.

 Then, it will go to the next stage. C & F change.


* Min. abnormal temp. alarm limit (0.0 ~ -5.0 °C Default Setting:-3.0 °C)





Then, it will go to the next stage, °C & °F change.

(The others, same with abnormal high temp, alarm function.)

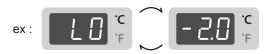

※ It shows that the set temperature is 2°C (2°F) lower than the
actual cabinet temperature of the machine.

* Just press (o.k) button if you do not need to change setting.

18




4. Function Settings

(3) How to change Centigrade & Fahrenheit / Work Automatic Pumping System by Force

How to change Centigrade & Fahrenheit

- ► From the abnormal LOw temperature alarm setting, press (o.k) button once and "C.F" (Celsius/Fahrenheit) will be displayed for 0.5 second and LED for "C" will blink.
- ▶ Select °C or °F with ◆ or → button.
- ► After selecting °C or °F , press (o.k) button to save. [Default setting : °C]



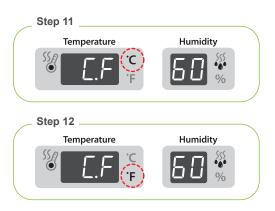

How to Work Automatic Pumping System(APS) by Force

- ▶ Press (+) button for about 10sec. then pump motor works for 2min. by force. It still works for 2min. even if you remove your finger from the button.



If you press button for about 10sec. pump will operate for about 2min. Press any key to remove.

5. Management

(1) How to Disassemble and Clean the Incubator

INSECT TO

How to Disassemble and Clean the Incubator





Be sure to take off power cord before cleaning the incubator.


► After incubation terminated, never leave the incubator covered with view window.

Moisture inside of the incubator vaporizes and it can stick to electric parts causing operation failure. Be sure to dry the incubator perfectly.


► Remove main controller from view window. [Refer to the page 9]

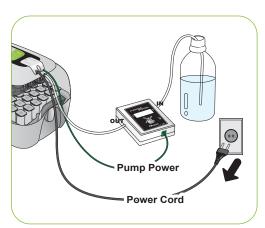

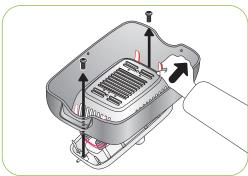

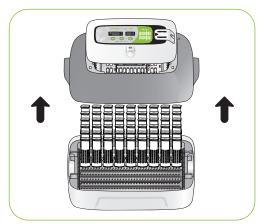

- Wipe the main controller smoothly with soft brush. Be careful not to give impact to sensor part or heating part.
- ▶ Temp. & Humidity sensor is assembled with connector for easy replacement in case of trouble.

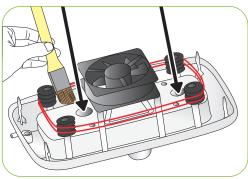


20

INSECT TO

5. Management

(1) How to Disassemble and Clean the Incubator

▶ Wash the window, tray, tray hatching pad, insect house and main body(bottom) with warm water.

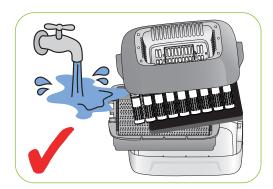


▶ Do not wash Automatic Pumping- System(APS) with water, just wipe with a soft cloth.



Do not use benzene or thinner when wipe the unit. It can cause transformation or decoloration.



▶ Wipe the exterior with a soft cloth, and completely dry before storing.

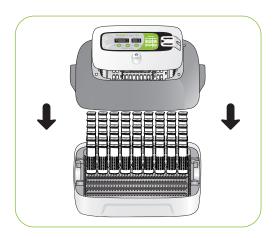


6. Product Information

(1) How to Replace the Fuse




How to Replace the Fuse


If power does not turn on even if you connect the plug, check fuse.

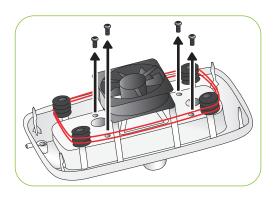

▶ Open the cover as Fig. then you can see fuse equipped PCB. Remove damaged fuse with screwdriver.

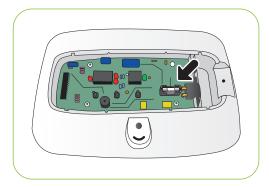

- ► Replace with new fuse. (250V 2.0A Ø5*20mm)
- ▶ Reassemble with the opposite order of disassemble.
- If the cover is not completely closed, moisture preventing ring will not attached correctly and moisture can be flow in to the PCB part.

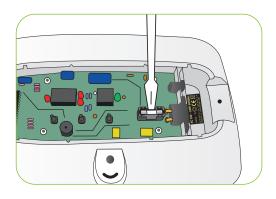

22



6. Product Information

(2) Troubleshooting

Troubleshooting

Circumstance	Items to be confirmed	Solutions		
No signs of power to the incubator.	 Check electric cord is properly connected. Check if there is a power outage. Check whether the plug is damaged. Check the fuse is not blown. [Refer to the page 22] 	 Connect the cord again. Check the main socket with any other electric appliances. Try to insert to another outlet. Replace the spare fuse in the fuse box. (250V 2.0A) 		
When temperature is too high / When temperature is too low Check your setting temperature. Check the air circulation FAN is working. Initialize the incubator. [Refer to the page 16]		 Set the temperature as required. Take the power plug out of the outlet, and disassemble the main body for cleaning around the air circulation fan with brush. 		
When humidity is too high / When humidity is too low	 ▶ Check water is enough. ▶ Check setting humidity. ▶ Initialize the incubator. [Refer to the page 16] <when low=""></when> ▶ Check if evaporating pad is fit well. ▶ Check if view window is closed completely. <when high=""></when> ▶ Check ambient conditions. ▶ Check if there is water in bottom body. 	 Supply water. Set required humidity. <when low=""></when> If evaporating pad does not fit well, humidification device will not work well. Close view window exactly. <when high=""></when> Adjust the conditions of room temperature [Refer to the page 8] Remover water inside of bottom body. 		
The machine makes some noise A little noisy occurs from the inside of circulation fan.	 Check whether you put anything on the incubator. Check if there is any vibration, or check the incubator set on a table. Check if there is any hatching debris such as feather or eggs-shell in the incubator FAN. 	 Ensure the incubator is not touching any other items and nothing is on top of the unit. Move the incubator to a flat and even surface. Unplug the power supply and disassemble the main body for cleaning the egg turning tray and gear. 		
Dew forms in the incubator. - Dew condensation is normal with high humidity incubation.	► Check if incubation room temperature is too low.	► Check conditions of incubation room, and adjust as required.		
When pumping is slow	► Check Silicon Tube	▶ Adjust APS speed hole with screw driver.		



Test the incubator to be sure that the incubator works without any problem before placing eggs in the incubator. Please ensure you remember how to use the machine and maybe re-read instructions.

6. Product Information

(3) Specification



Specification

1-1 Name: Rcom Insect MAX 70

1-2 Model: MX-N70

Power AC 100[V]~120[V], 50/60[Hz] / AC 220[V]~240[V], 50/60[Hz]		
Power Consumptio	n Average 25[W] Max. 48[W]	
Temp. limits	20 ~ 42[°C] / 68 ~ 107.6[°F]	
Humidity Limits	20[%] ~ 70[%] (It can be changed according to the conditions)	
Weight	2.4[Kg]	
Size	(W)370×(L)231×(H)209[mm]	
Fuse Standard	250[V] 2.0[A] (Ø5*20[mm])	

You will be asking for service charge if you are in below cases.

- 1. Broken or Problem caused by 'No cleaning' or 'Improper handling' 2. Broken or problem caused by 'Not using distilled water'
- 3. Broken or problem caused by 'Trying Bird hatching in incubator' (Please don't try bird hatching in incubator. Please use Hatcher during bird hatching.)
- 4. Broken or problem caused by 'Wrong controling(Using)'5. Broken or Problem caused by using the machine out of User's Manual.
- 6. Broken or problem caused by 'Customer's mistake or fault'
- * Register your product Serial No. on our website for free 2 years gurantee.
- * Refer to website for how to register. (Free 1 years gurantee in case not registerd.)
- Product Registration Procedures

If you are a new member of Rcom, you need to log in our website

- 1. Click "SIGN UP" on top of the right corner.
- 2. Fill out the information on the blanks.

If you already registered your information on the Rcom website, please log in

- 1. Click "Login" on top of the right corner.
- 2. Select "Customer" and click "Register Products".
- 3. Fill out the information on the blanks.

You will need below information to register on our website.

- 1. Your name
- 2. Your (home/company) address & Email address, country
- 3. Your product(model) name & Serial number
- 4. Date of purchase & Name of purchased shop you bought from

24

North America Sales & Service www.RcomCo.com

Rcom Incubator Series



Rcom PRO MINI (PX-03)



Rcom KINGSURO MAX 20 (MX-SURO)



Rcom MAX 20 (MX-20)



Rcom PRO 50 (PX-50)



Rcom REPTILE MAX 90 (MX-R90)




Rcom REPTILE PRO 90 (PX-R90)



Rcom REPTILE MAX 230 (MX-R230)



Rcom REPTILE MAX 230 (PX-R230)

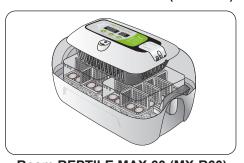

Rcom REPTILE MAX 60 (MX-R60)



Rcom DIGILOG (PX-V1)



Rcom DIGILOG2 (PX-V2)

North America Sales & Service www.RcomCo.com

Rcom is designed for user's easy and convenient use. It can be changed without notice for improvement in performance, design, treatment, and software, etc.



www.RcomCo.com



North America Sales & Service

www.RcomCo.com

MANUFACTURER

AUTOELEX CO.,LTD.

• OFFICE: 218, Sangdong-ro, Sangdong-myeon, Gimhae-si, Gyeongsangnam-do, 50805 Korea

Rcom is the Best Design with High Quality, Value, and Performance!

Rcom Insect Incubator Max 70 User Manual